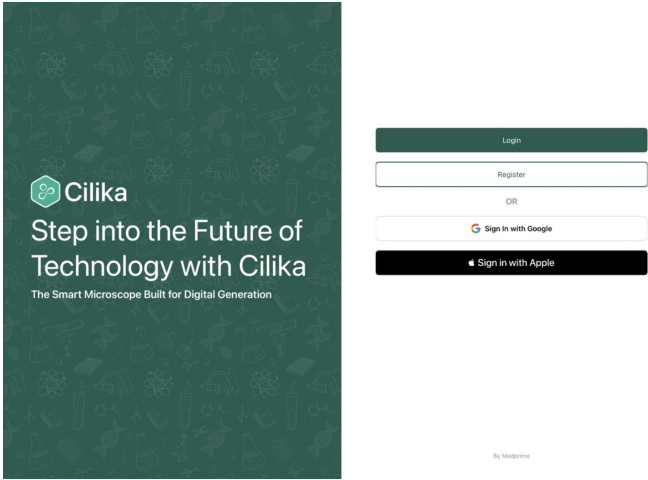
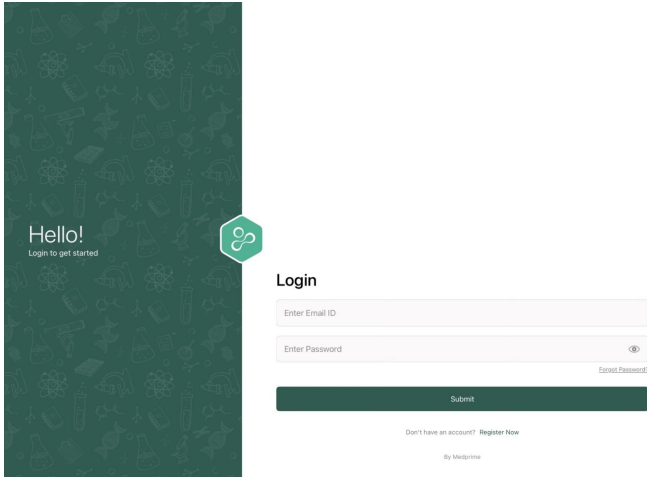
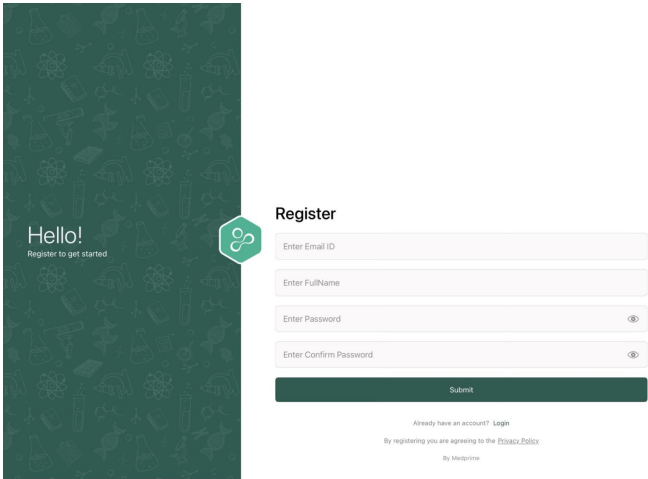


Initial Configuration

1. Install the "Cilika Reticulocyte" application from the "App Store"
2. Once the application is installed, the user can find the application in the home screen at the bottom.
3. The application starts with **Register/Login Screen**.
4. Login is for the users who have already used *cilika* in the past and have an account.
5. Registration is for the users who haven't used *cilika* in the past.
6. Registration requires the user to enter their email, their full name, and a new password.

 The login screen features a dark green background with a pattern of scientific icons. On the left, the Cilika logo is displayed above the text "Step into the Future of Technology with Cilika" and "The Smart Microscope Built for Digital Generation". On the right, there are four buttons: "Login", "Register", "OR", "Sign in with Google", and "Sign in with Apple". At the bottom, it says "By Medprime".	 The login screen features a dark green background with a pattern of scientific icons. On the left, the text "Hello! Login to get started" is displayed next to the Cilika logo. On the right, there is a "Login" section with two input fields: "Enter Email ID" and "Enter Password". Below these fields is a "Submit" button. At the bottom, it says "Don't have an account? Register Now" and "By Medprime".
 The register screen features a dark green background with a pattern of scientific icons. On the left, the text "Hello! Register to get started" is displayed next to the Cilika logo. On the right, there is a "Register" section with four input fields: "Enter Email ID", "Enter FullName", "Enter Password", and "Enter Confirm Password". Below these fields is a "Submit" button. At the bottom, it says "Already have an account? Login" and "By registering you are agreeing to the Privacy Policy" and "By Medprime".	

Revision #5

Created 2026-07-21 16:53:45 IST by Rashi

Updated 2026-07-28 15:36:02 IST by Rashi