

App Settings

1.5 App Settings

Go to the back to the home page, and tap on the settings icon. A pop up window appears on the screen with various settings option.

Image Compression

1. **This setting offers the user to select, image quality type.**
2. **Low: Medium: Original:**

Video Compression

1. **This setting offers the user to select, image quality type.**

Screen Recording

Allows users to record live view screen, can be used in demonstrations and trainings.

Horizontal Flip

1. **Tap on the Horizontal Flip toggle to flip the live view screen image horizontally.**

Vertical Flip

1. **Tap on the Vertical Flip toggle to flip the live view screen image vertically.**

Measurement Unit

1. **Tap on the toggle to change the measurement unit to either micrometer or millimeter**

Enable Tiff Format

1. **Tap on the Enable Tiff to stores the image in the highest quality.**

Lens Profile

1. **Go to lens profile by tapping on it**
2. **A default lens profile will already be active.**
3. **Scroll down to set up another lens profile.**
4. **Add your custom lens and give it a personal color.**
5. **Tap on the plus icon on the write to add it.**

6. **Tap on the lens you want to see on the live screen. Selected lenses will turn black. Preview the lens you have selected.**
7. **Add a profile name and tap on add. Lens profile is added.**
8. **Scroll back up and tap on your profile to activate it.**
9. **To delete the lens profile, tap on the edit option above the status.**
- 10.

Object Storage

Quick Share

1. **Helps direct file transfers between your iPad and a desktop computer (Windows or Mac) within the same local network.**

Reference Image

1. **An external reference image can be added on the live view screen and can be positioned according to the users preference.**

Watermark General

1. **Watermark can be used as an identifier.**
2. **The watermark can be customized according to the**

Watermark Cilika

Zoom

Scale Setting

Auto Hide

RBC Template

hide Calibration

Enable Grid

Unit

Size

Opacity

Line Type

Color

Show frame

Revision #9

Created 2026-05-20 13:23:59 IST by Rashi

Updated 2026-05-23 15:34:37 IST by Rashi